



Colifast FAST – Quick guide

Equipment needed:


FAST Vial

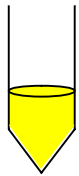

Sterile Scalpel

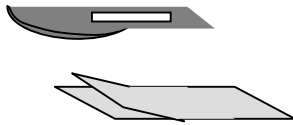


Incubator

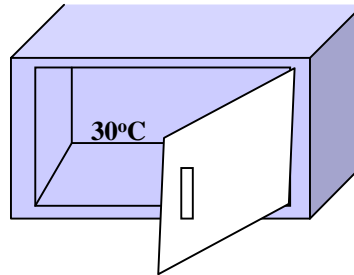

Clock/Timer



Procedure:

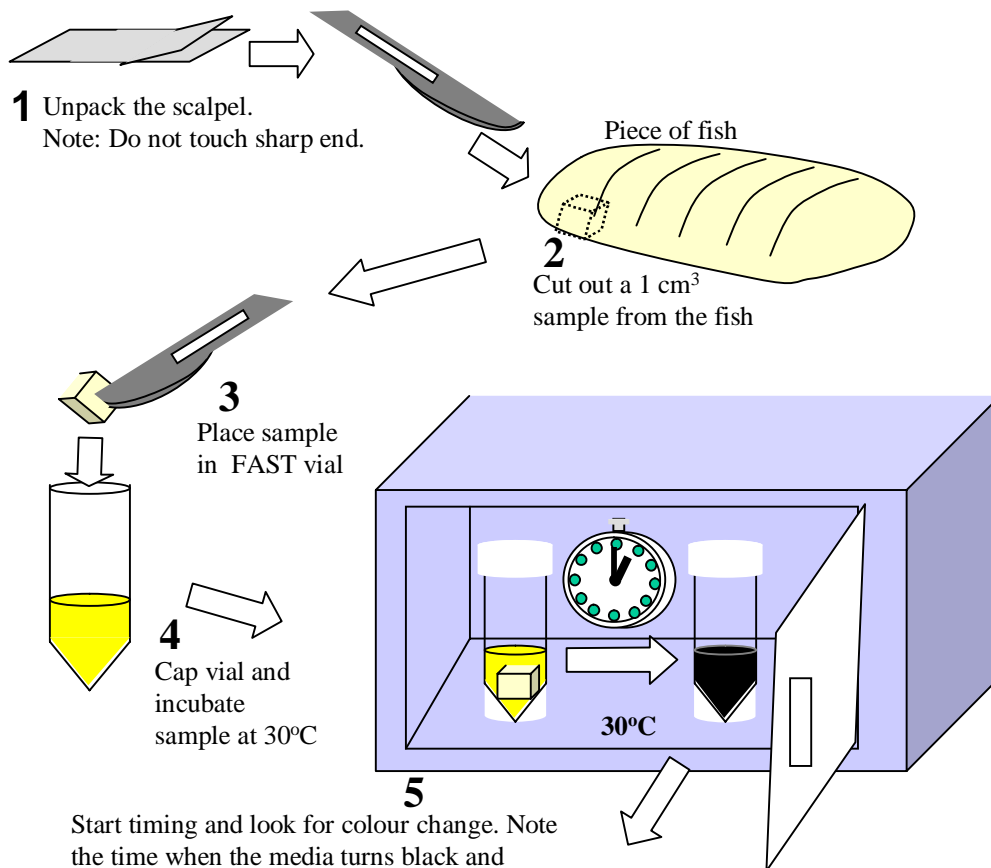

Note: Beware contamination during sampling. Do not touch the blade of scalpel or sample. Keep vials sealed. Use scalpel only once.

Hours to colour change Yellow → Black	Sulphide Producing Bacteria/g	Quality Guideline
3-6	>5 000 000	Bad
7-10	500 000 - 5 000 000	Marginal
11-12	1000 - 500 000	Good
>12	<1000	Very good

COLIFAST®